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Revision changes

06-05-2019 Time until preparation is changed in time until freezing.

06-05-2019 REMP-tube is change in MATRIX tube.

05-07-2021:

- SOP Format is changed.
- Pre analysis conditions are added.
- Concentration in ng/ul instead of ug/ml.
- Specifications Matrix tube added.
- Removed: After 5 issuances, make a new working solution from the stock solution at the time that the sixth issuance is requested.
- Origin is added to data management.
- Making of working solution the first time the DNA is issued is removed from this SOP and is never done.
- Added: If DNA concentration less than 30 ng/ul the isolation is failed, see Deviations.

1. Goal

To describe how DNA from saliva must be prepared and stored for purposes of the Radboud Biobank (RB), so that:

1. Interested parties including potential users know how the biomaterial has been handled.
2. The laboratory can assess whether they can process the biomaterial as described under paragraph 2 "Preparation".
3. The laboratory knows the requirements for the collection, processing and storage of the biomaterial to enable registration of deviations.
4. The sub-biobank or project knows the requirements for the collection and transport of the biomaterial to ensure correct delivery.
5. The RB is able to couple this procedure to the biomaterial in storage to determine the fitness for purpose.

1.1 Scope of application

This procedure is applicable to all employees concerned at the RB, the Sub-biobank/Project and the Department of Human Genetics of the Radboudumc.

2. Protocol isolation and storage of DNA from saliva

Collection

Preanalysis

Conditions for the collection of saliva for DNA investigations.

- No eating, drinking, smoking, cleaning of teeth or use of mouth wash or chewing gum 30 minutes before collection of saliva.
- Remove lipstick if present.
- Before collecting, rinse mouth out with water and then wait 5 minutes.

Delivery

In collection tube.

Collection

2 ml saliva (or max. 5 "sponges" from Oragene) added to Oragene/saliva vial (Oragene kit; Oragene DNA (OG-500) from DNAgenotek, Ontario, Canada).

Temperature

Keep tubes at room temperature until isolation.

Preparation

Time until freezing

Isolate the DNA as soon as possible.

Title: SOP Isolation and storage of DNA from saliva

One year maximum (shelf-life of the saliva sample after combining with the Oragene-solution).

Isolation protocol DNA isolation from Saliva with ChemagicStar ([Qdoc 084294](#)).
 Estimated yield of approximately 80 µg DNA.

DNA-concentration and volume After isolation determine the DNA-concentration (in ng/ul) and the OD-ratio 260-280 nm from the stock solution and register in the Biobank management system.

Storage

Aliquotation Store the DNA in a 500 ul stock solution (1 aliquot).

Type of tube MATRIX-tube (Matrix™ 2D Barcoded Open-Top Storage Tubes, ThermoFisher Scientific).

Temperature -20°C (Liconic).

Miscellaneous

Data management Register the following data in the biobank management system:
 1. Date the saliva was collected.
 2. Input-date is time of storage at -20°C after DNA-isolation.
 3. Identification number (pseudomised).
 4. Name of sub-biobank/ pearl/ project.
 5. Fraction number.
 6. OD-ratio and DNA-concentration.
 7. Origin of the sample.

Deviations Note all deviations from this procedure in the Biobank management system. If the DNA concentration is less than 30 ng/ul the isolation is failed. Failed isolations must be registered, and the RB must be informed. The stocks are stored in the Liconic.

3. Quality assurance

Review takes place via i) an internal audit (see the [audit year plan](#) in DMS) and ii) 2-monthly performance reviews of deviations registered in the biobank management system (DMS map "[Registraties en controles](#)").

3.1 Performance indicator

95% of the sub-biobank samples have been prepared and stored within the parameters of this procedure.

4. Comments

Although this SOP is used for collection of Parelsnoer DNA, this SOP differs from the standard of Parelsnoer. Parelsnoer wants a minimum of 2 and prefers 3 or 4 aliquots. Given the costs and the frequency DNA is issued the RB decided to store only 1 DNA stock per isolation.

5. Relevant documents

Qdocs, literatuur, wet en regelgeving, website e.d.

Qdocs, literature, legislation and codes of conduct, website etc.

link	Title
Qdoc 083613	Monstermanagement SOP KGCN

Qdoc 045505	PVS 10.05 Primair laboratoriumproces, DNA-isolatiefaciliteit
Qdoc 072370	DNA_NORM_ALLIN
Qdoc 084294	DNA_isolatie_ALLIN
Qdoc 070512	<u>Liconic Bio.Li.X</u>
Qdoc 083792	Liconic vriesrobot, werking koelsysteem en alarmering
Qdoc 069283	Bewaartermijnen Genetica

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